



The SMA OneTouch[®] Integrated Control System (ICS) is a computerized, automated viable air monitoring system that controls calibrated, precise air sampling through individual or multiple SMA Atriums[®]. SMA Atriums are stainless steel collection devices that direct air from the environment to impact onto a media plate. The media plate is then evaluated to determine the amount of viable contamination in the air sample.

The SMA OneTouch ICS strictly regulates the air flow rate throughout a sample cycle. The system automatically alarms the operator visually and audibly if the sample becomes compromised or is aborted. In addition, the SMA OneTouch ICS continuously monitors the air flow rate while sampling and will alert the operator if the flow rate deviates from the established 1 CFM flow rate.

The SMA OneTouch ICS has a touchscreen interface that incorporates facility floor plans and SMA Atrium locations for sample monitoring. User accounts have access levels which limit their access to specific functionality.

VAI also offers the SMA OneTouch ICS for Isolators. This SMA OneTouch ICS configuration includes all the aforementioned benefits and provides secure sampling inside the isolator environment.

Product Uses

- Program, start, abort, and monitor all sampling through integrated SMA Atriums
- Monitors all sample parameters (e.g. sample volume, sample duration, and vacuum pump status)
- Provides visual indication of all activity throughout the sampling process
- Immediate audible and visual alarms will sound if the sample is compromised or aborted

Features and Benefits

- Installation requires minimal tubing and wiring which simplifies installation into new and existing facilities
- Can be integrated with existing facility monitoring and data collection systems using industry standard networking
- Provides system-wide viable monitoring capabilities from a single interface
- Can be remotely monitored using a variety of handheld, tablet, and computer-based devices

Proven Solutions for Viable Air Monitoring

SMA OneTouch ICS Interface

Description	The Interface is a touch screen that is used to monitor and control the SMA OneTouch ICS and all integrated SMA Atriums.
Enclosure(s)	316L or 304 stainless steel (depending on configuration)
Input power	24 VDC power sourced from the Controller
Size	Interfaces are available in multiple sizes ranging from 6.5 to 15 inches.

SMA OneTouch ICS Controller

Description	The Controller is a computerized system which incorporates a Programmable Logic Controller (PLC) and powers the Interface and Flow Center. It automates all system monitoring, sampling, and alarming functions.
Enclosure(s)	316L or 304 stainless steel, free standing
Input power	120 VAC (North America), dependent on system design

SMA OneTouch ICS Flow Center

Description	The Flow Center contains a series of high accuracy mass-flow controllers (1.0 CFM \pm 2%) that automatically regulate the air flow on each sampling line.
Enclosure	316L or 304 stainless steel (depending on configuration)
Input power	24 VDC power sourced from the Controller

Tubing and Cable Connections

Cables	Category 5e networking cable (Cat 5e) or higher
Tubing	¼ and ⅜ inch ID Hytrel lined PVC tubing ⅜ inch ID Vardex reinforced vinyl tubing

SMA Atriums

Description	An SMA Atrium is a stainless steel device that directs air from the environment to impact onto a media plate. The media plate is then tested to determine the amount of viable contamination in the environment.
-------------	--