


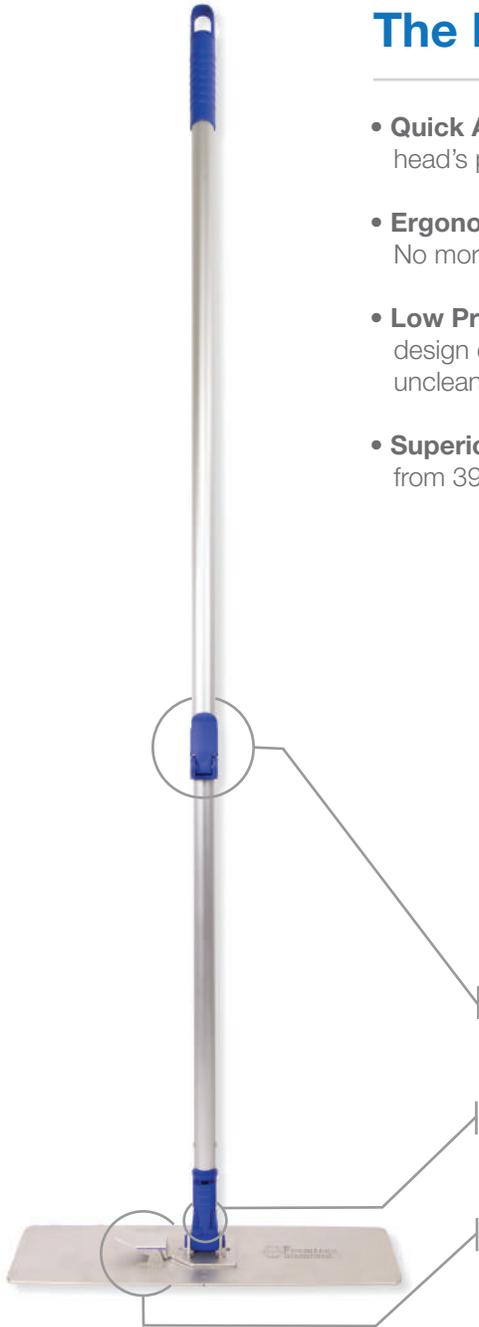

## The PharmaMOP® GO Difference

- **Quick Assembly & Disassembly.** Head frame easily slides into the mop head's pockets for fast insertion and disposal of mopheads.
- **Ergonomic Foot-Switch.** Discard mop heads while standing upright. No more touching soiled mops while discarding — No more stooping.
- **Low Profile Design.** A two-direction 360° swivel with a 1.5" low profile design enables unsurpassed access to tight spaces reducing previously uncleanable "dead spots".
- **Superior Handle.** Fully autoclavable, lightweight, aluminum handle extends from 39-65" (99-165cm) making it ideal for use on walls, floors & ceilings.



## Features & Benefits

- 1** Extendable Handle provides efficient cleaning of walls, floors and ceilings.
- 2** 360° Swivel Lock allows for both Figure 8 and Flat Mopping.
- 3** V-Lock Lever provides easy, quick insertion & disposal of mop heads.



Foamtec International  
720 Venture Dr.  
Waco, TX 76712

Foamtec International WCC Division  
1-760-599-7160

www.foamtecintlwcc.com | customerservice@foamtecintlwcc.com



## Sahara+ PharmaMOP® GO Mop Head



- **Advanced Construction.** An absorbent cleanroom foam core is ultrasonically welded between a layer of a non-linting hybrid microfiber and laminated with dual Sahara® Foam Scrubbing Strips for exceptional cleaning.
- **Minimize Surface Microbial Contamination.** Unique mop head materials dislodge, entrap and remove residues and soils while evenly applying disinfectant to meet contact time requirements.
- **Reduce Work Time.** Ultrasonically sealed pockets allow the mop frame to slide into the mop head and be ready-for-use within seconds.
- **Ultimate Compatibility.** Provides efficient cleaning of walls, floors and ceilings. Compatible with bleach, QUAT, peracetic acid, peroxide, and IPA-based cleaning and disinfecting agents.
- **Quality Manufacturing.** Manufactured in an ISO 13485 & ISO 9001 facility on a fully automated fabrication line. Ideal for use in ISO 3-9 cleanrooms. Available gamma irradiated, SAL 10<sup>-6</sup>.

## Features & Benefits

- 1** Sahara® Foam Scrubbing Strips dislodge contaminants and provides superior abrasion resistance.
- 2** Lint free hybrid microfiber fabric entraps & removes microbial contamination and stubbornly adhered disinfectant residues.
- 3** The combination of Sahara® Foam Scrubbing Strips and Microfiber provide optimal gliding even on high friction, non-skid cleanroom floors.
- 4** Encased Foam core provides high absorbency to evenly apply disinfectants & 70% IPA to enhance cleaning performance and to meet contact times.
- 5** Ultrasonically Sealed Edge provides unsurpassed in-use particulate control.



Foamtec International  
720 Venture Dr.  
Waco, TX 76712

Foamtec International WCC Division  
1-760-599-7160

www.foamtecintlwcc.com | customerservice@foamtecintlwcc.com



PharmaMOP® GO

# Bucket & V-Bracket

*Hands-free, Fiber-free Cleanroom Mopping*

## PharmaMOP® GO Bucket & V-Bracket

The stainless steel v-bracket inserts into a specially-designed mop bucket enabling a hands-free assembly of the mop head onto the mop frame. In addition, mop heads can be pre-saturated in the bucket for cleaning SOP's that require single-use mop heads.

The bucket and bracket system accommodate up to 12 mop heads. All components are autoclave safe and are designed to be used in ISO Class 3-9 cleanrooms.



Foamtec International  
720 Venture Dr.  
Waco, TX 76712

Foamtec International WCC Division

1-760-599-7160

[www.foamtecintlwcc.com](http://www.foamtecintlwcc.com) | [customerservice@foamtecintlwcc.com](mailto:customerservice@foamtecintlwcc.com)

## PharmaMOP® GO Technical Data

### Recommended Autoclave Settings\*

Exposure Temperature	121°C
Exposure Time	30 Minutes
Cycle Type	Gravity
Cycle Time	20 Minutes
Pressure	17.5 psig

PharmaMOP® GO Mop Head	Particle Data via Helmke Drum Testing (Particles/Mop Head)
0.5 - 10.0µm	≤ 700

\* The recommended autoclave settings are intended for the Lever Lock Mop Handle, Head Frame & Cleanroom Mop Heads.

## PharmaMOP® GO Ordering Information

### PharmaMOP® GO Mop Heads

Part No.	Description	Mop Heads/Pack	Packs/Case
FS879SH-1	Sahara+ PharmaMOP® GO Mop Head - Individually Packed	1	200
FS879SH	Sahara+ PharmaMOP® GO Mop Head	5	40
FS879SHST	Sterile Sahara+ PharmaMOP® GO Mop Head - Individually Packed	1	200

### PharmaMOP® GO Handles & Accessories

Part No.	Description	Quantity
FT430	15" PharmaMOP® GO Handle w/ Locking Head Frame with VLock on Telescoping Handle	1
FT432	Stainless Steel V-Bracket for Hands Free Assembly and Disassembly (15.8" x 5.5" x 7.7")	1
FT408-DB	Double Bucket System with Frame	1
FT409BL	Blue Cleanroom Bucket (20.5" x 10.5" x 10.5")	1

### PharmaMOP® GO Starter Kit

Part No.	Description	Quantity
FS4500-PHMGO	PharmaMOP GO Starter Kit - FT430 (1), FT432 (1), FT409BL (1), FS879SH (25)	1



**Foamtec International**  
720 Venture Dr.  
Waco, TX 76712

**Foamtec International WCC Division**

**1-760-599-7160**

www.foamtecintlwcc.com | customerservice@foamtecintlwcc.com